

# COLORIMETRIC PH TEST EQUIPMENT


#### **CODE 5858**


#### **USE OF THE OCTET COMPARATOR**



The Octet Comparator contains eight permanent color standards. A test sample is inserted into the openings in the top of the comparator. The sample can then be compared to four color standards at once, and the value read off the comparator. For optimum color comparison, the comparator should be positioned between the operator and a light source, so that the light enters through the special light-diffusing screen in the back of the comparator. Avoid viewing the comparator against direct sunlight or an irregularly lighted background.

## **pH TEST PROCEDURE**

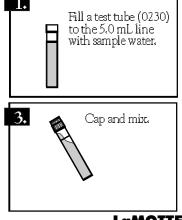

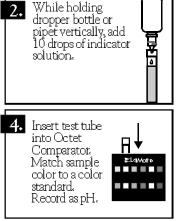

## LaMOTTE COMPANY

## Helping People Solve An alytical Challenges®

PO Box 329 • Chestertown • Maryland • 21620 • USA 800-344-3100 • 410-778-3100 (Outside U.S.A.) • Fax: 410-778-6394 Visit us on the web at www.lamotte.com